


Rotavapor® R-250 EX Technical data sheet

The Rotavapor R-250 EX is the EX rated version of the Rotavapor R-250. It fulfills the high requirements of the ATEX regulations and fits perfect to your Ex-proof production laboratory. It is designed for easy and safe operation combined with high performance.

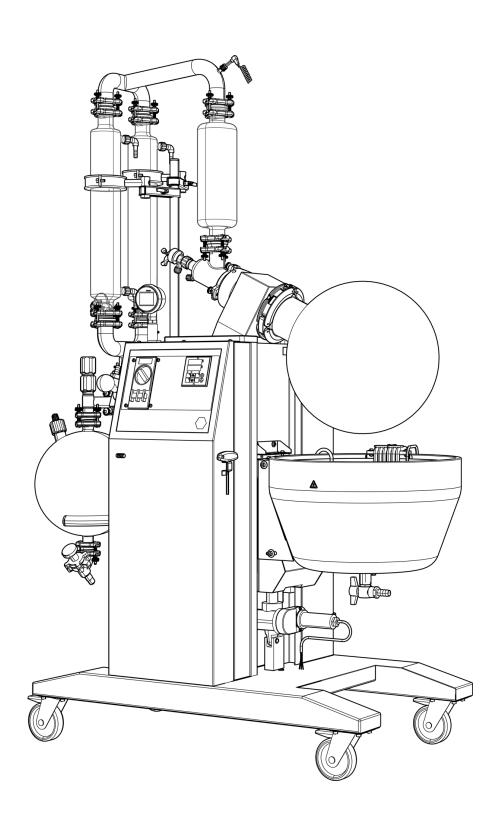


Scope of delivery

All Rotavapor R-250 EX models are delivered ready to use, including all needed tools and are complete of:

Components	R-250 EX
Vacuum controller UDS7 EX	1
Vacuum valve EX	1
Glass configuration according order	1
Evaporating flask according order	1
Vacuum hose to the vacuum source (PVC), 3.5 m	1
Vacuum hose to the vacuum controller (rubber), 3 m	1
Cooling water hose 19/26 (PVC), 2 m	1
Cooling water hose 14/18 (PVC), 3 m	1
Cooling water hose 10/14 (PVC), 2.2 m	1
Bath replenishment hose (PVC), 2 m	1
Feed hose (PTFE), 2 m	1
Seal removal tool	1
Operation manual	1

Order code

Choose the configuration according to your needs:

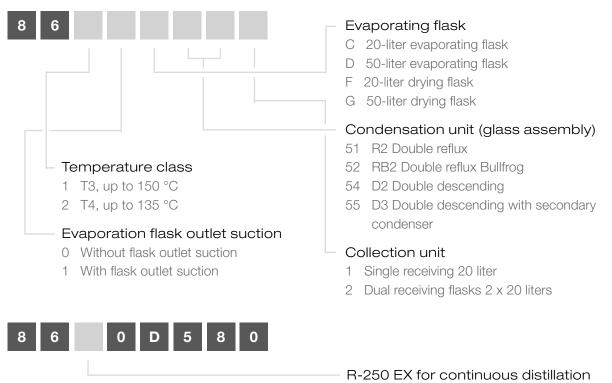

- 1 Version T3 up to 150 °C
- 2 Version T4 up to 135 °C

Dimensions and weight (hight) *226 cm *210 cm

D2	D3	RB2	R2
Low boiling points and/or foaming products		High boiling points	
	Minimum emissions	Reflux	x reactions

Reduced height

	Dimensions (W x H x D)	Weight
Assembly R2	1410 x 2300 x 830 mm	285 kg
Assembly RB2	1410 x 2100 x 830 mm	285 kg
Assembly D2	1410 x 2260 x 830 mm	290 kg
Assembly D3	1410 x 2260 x 830 mm	295 kg

Technical data

Electrical

Voltage	400 – 440 V (3Ph, N, G)
Frequency	50 – 60 Hz
Max. power consumption	7500 – 9100 W

Standards

EX label	Ex de ia [ia] IIC T3 and T4
CE marking	Yes

Safety

Salety	
Safety coated glassware	No, due to the risc of electrostatic charges
Over temperature protection of the bath	Separate monitoring circuit with manual reset Error if temperature is 15 °C above set value
Rotation	Soft start
At any Error	Heater off, rotation off, bath lowers Indication of error

Display

Bath temperature	Yes (1 °C steps)
Vapor temperature	Yes (1 °C steps)
Set rotation speed	Yes (1 rpm steps)
Set bath temperature	Yes (1 °C steps)
Actual vacuum	Yes (1 mbar steps)
Set vacuum	Yes (1 mbar steps)

Cooling

Consumption	200 – 400 liter/hour (adjustable via integrated needle valve)
Restriction	max. 2,7 bar abs. without pulsation

Heating (bath)

Medium	Water or oil
Temperature range	Ambient to 115 °C (T4) or 150 °C (T3)
Heating capacity	6600 W (3 W/cm²)
Accuracy	+/- 2 °C

Rotation

Motor	400 V (3Ph; 2A at 50 Hz)
Controlling	Frequency inverter
Speed	0 – 100 rpm
Accuracy	+/- 1 rpm at 5 rpm up tp +/- 5 rpm at 100 rpm


Materials

Gear headAluminum cast (3.2373)PaintingPowder coated with Epoxy (EPX)Bath panStainless steel 1.4404 (AISI 316L)Heating elementStainless steel 1.4404 (AISI 316L)GlassBorosilicate 3.3In contact with productFDA compliant materialsLeakage of the complete system< 1 mbar/min	Housing	Stainless steel 1.4301 (AISI 304)
Bath pan Stainless steel 1.4404 (AISI 316L) Heating element Stainless steel 1.4404 (AISI 316L) Glass Borosilicate 3.3 In contact with product FDA compliant materials	Gear head	Aluminum cast (3.2373)
Heating element Stainless steel 1.4404 (AISI 316L) Glass Borosilicate 3.3 In contact with product FDA compliant materials	Painting	Powder coated with Epoxy (EPX)
Glass Borosilicate 3.3 In contact with product FDA compliant materials	Bath pan	Stainless steel 1.4404 (AISI 316L)
In contact with product FDA compliant materials	Heating element	Stainless steel 1.4404 (AISI 316L)
	Glass	Borosilicate 3.3
Leakage of the complete system < 1 mbar/min	In contact with product	FDA compliant materials
	Leakage of the complete system	< 1 mbar/min

Sensors

Vapor temperature	PT-1000, 2 wire
Bath temperature	PT-1000, 2 wire
Vacuum	Piezoresistive measuring cell; stainless steel 1.4435 (AISI 316L)

Performance

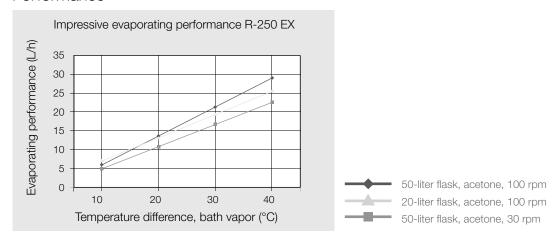

For other solvents the following maximum general figures apply:

Water	Methanol	Ethanol	Toluene	Ethylacetate	Hexane	Trichlorethane
5.8 L/h	14.0 L/h	19.5 L/h	36.5 L/h	38.0 L/h	55.0 L/h	40.5 L/h

Accessories

	Order number
Vacuum pump MD4C EX PTFE membrane pump with a capacity of 3.7m3/h and a final vacuum of 3mbar. 230 V / 50 Hz	040197
Manual flask handler 50 liter For easy mounting and removal of the 50 liter evaporating flasks along with safe transport.	041414
Manual flask handler 20 liter For easy mounting and removal of the 20 liter evaporating flasks along with safe transport.	041410
Flask crane (EX version) For the safely secured transport of a 50 liter flask. Incl. the 50 liter manual flask handler	041493
Vapor duct with sinter plate With frit P3 protects the condenser assembly against powder and dust during the drying process.	041100
IQ / OQ	40786